

Status Bar Icons: Internal vs. Network Storage

PATsmart™ REBEL® System

Support Document

Below are the status bar icons that are listed in the PATsmart™ REBEL® System interface and their associated meanings. For more information, visit the Repligen [Knowledge Base](#) or contact TechSupport@repligen.com



Internal Storage Configured (REBEL is ready to run in this mode.)



Network Storage Configured (REBEL is ready to run in this mode.)



Files Syncing (REBEL has files stored locally that are transferring to the network.)



Network Locations Unreachable (REBEL is configured for network storage, but the read/write directory is not accessible.)

- To resolve, navigate to read/write directory > Test Permissions > Apply.

Files Queued for Export (REBEL has files stored internally that will be sent to the network when the connection is reachable before a new batch begins.)

- To resolve, navigate to read/write directory > Test Permissions > Apply.

Customer Service

Repligen Corporation
685 Route 202/206
Bridgewater, NJ 08807

TechSupport@repligen.com

(908) 707-1009

repligen.com

© 2025 Repligen Corporation. All rights reserved. The trademarks mentioned herein are the property of Repligen Corporation and/or its affiliate(s) or their respective owners.

