

Raman-Based Lactate Feeding Impact on Culture Characteristics and Product Quality in Mammalian Cell Cultures

Stephanie R. Klaubert¹ • Awab Nehela¹ • Stephen P. Driscoll² • Morgan Siegmann² • Christopher D. Brown² • Scott E. Miller²
¹908 Devices Inc., Boston, MA, USA, ²Repligen, Marlborough, MA, USA



Repligen Corporation now owns the life sciences PAT product portfolio of 908 Devices Inc. Please contact Repligen for further inquiries.

Motivation

The biopharmaceutical industry is increasingly motivated to improve process control, understanding, and repeatability using process analytical technologies (PAT). More specifically, Raman-based optical, in-line measurement methods have shown promise as a way to monitor critical performance parameters, including glucose and lactate concentrations, in real-time. Recently, this has translated to the implementation of Raman-based glucose feeding strategies to modulate productivity and product quality in Chinese hamster ovary (CHO) cell cultures.¹ The current work implements a novel Raman-based solution that does not require any chemometric models for glucose, lactate and biomass measurements. The PATsmart™ MAVERICK® System was directly, without any training in the specific bioprocess, deployed to **dynamically control lactate feeding**. We compared feeding strategies as means of modulating product quality and improving culture conditions, primarily through decreased ammonia accumulation.

Methods

Experiment Conditions

Bioreactor

Distek BIOne 1250 3 L; Biostat A plus 3 L

Cell Line

NIST-CHO mAb

Seeding Density

0.5x10⁶ cells/mL; 4.5x10⁶ cells/mL

Medium

Ex-Cell Advanced CHO Fed-batch Medium (SAFC)

Feed

Ex-Cell Advanced CHO Feed 1 (SAFC) w/o glucose fed at 5% EOD

Glucose Feeding Strategy

Bolus feed with threshold <4 g/L, supplement to 6 g/L

Lactate Feeding Strategy

- 1.5 g/L lactate set point after consumption begins
- No lactate feed

Analytics

- In-line:** MAVERICK Raman-based real-time glucose, lactate, and biomass
- At-line:** Vi-CELL (Beckman Coulter), BioProfile FLEX2 (Nova Biomedical), and REBEL for amino acids
- Off-line protein quality attributes:** ZipChip coupled to Orbitrap Exploris 240 Mass Spectrometer (Thermo Scientific)

Experiment Goals

Goal

- Raman-based measurements used for dynamic lactate feeding for decreased ammonia accumulation in CHO cell culture
- Comparing **1.5 g/L lactate set point** to **no lactate feed**
- Both conditions had glucose bolus feed



PATsmart™ MAVERICK®
In-line glucose monitoring and lactate control




PATsmart™ REBEL® XT
At-line amino acids




PATsmart™ ZipChip®
CQAs

Growth and Productivity

Viable Cell Density

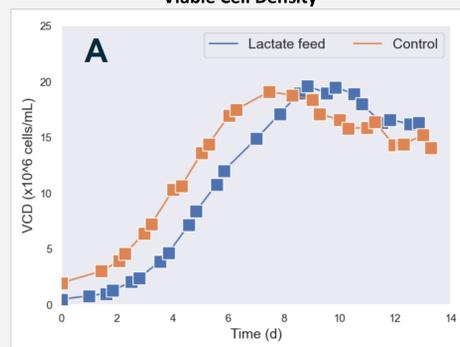

Titer

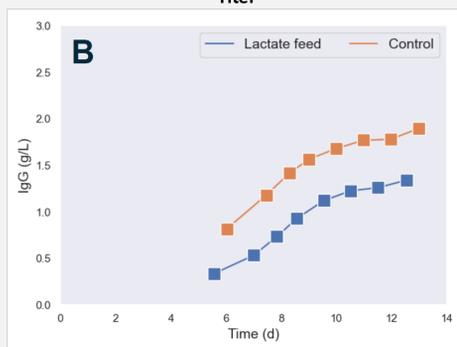

Cell Specific Productivity


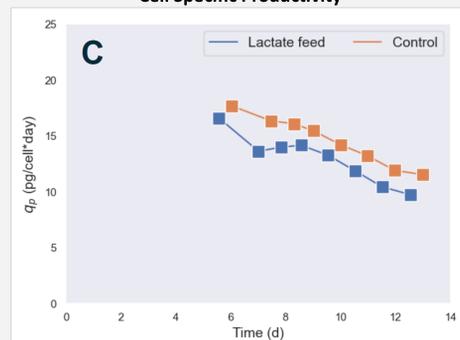



Figure 1. Lactate feeding impact on growth and productivity: A) Viable cell density (VCD), B) Titer, and C) Cell specific productivity (q_p) for control (orange) and lactate fed (blue) conditions. The lactate feed resulted in **comparable max VCD**, but **lower final titer and q_p** likely due to lower seeding density of the culture.

Metabolites

Glucose

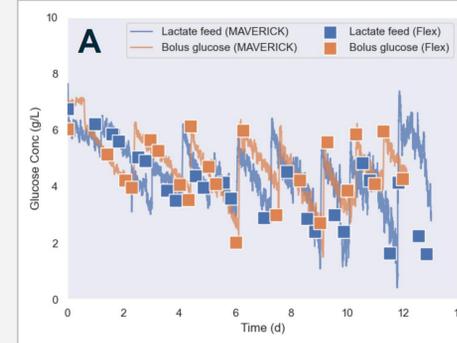

Lactate

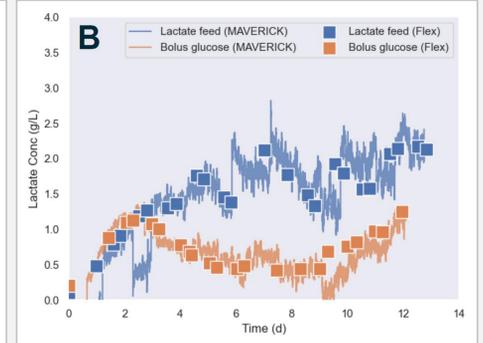

Ammonium

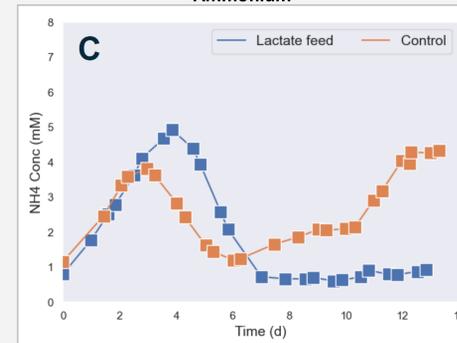


Figure 2. Lactate feeding impact on metabolites: A) Glucose, B) Lactate, and C) Ammonium concentrations for control (orange) and lactate fed (blue) conditions. The **lactate feed successfully reduced the ammonium accumulation** at the end of the culture.

Product Quality

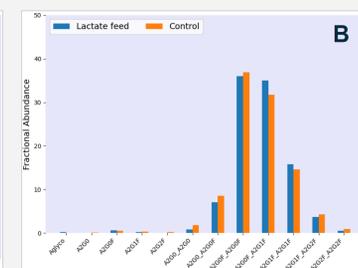
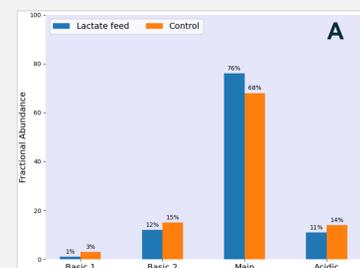


Figure 3. Lactate feeding impact on product quality: A) Charge variants and B) glycosylation for control (orange) and lactate fed (blue) conditions. The **lactate feed resulted in a higher percentage of mAbs in the main peak in the charge variant profile**, as well as a **higher abundance of more mature (G1 and G2) glycoforms**.

Alanine Profiles

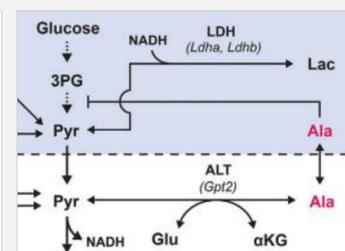
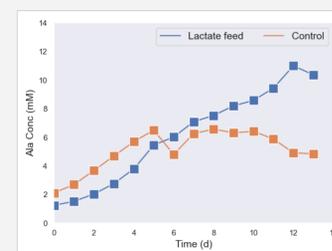


Figure 4. Lactate feeding impact on alanine levels: Sufficient lactate supply in the lactate fed cultures blocks a switch to alanine consumption. Instead of metabolizing alanine, lactate is used as a pyruvate source, leading to reduced levels of free ammonium from alanine metabolism.

Conclusions

- Easy and instant implementation of lactate control with MAVERICK
- Improved cell culture performance: ammonium profile
- Improved protein quality attributes: more mature, heterogeneous quality attribute profiles

References

- Domján J, Fricska A, Madarász L, et al. Raman-based dynamic feeding strategies using real-time glucose concentration monitoring system during adalimumab producing CHO cell cultivation. *Biotechnology Progress*. 2020.



Learn more at repligen.com