

# A Novel Integrated Strategy for the Detection and Quantification of the Neurotoxin $\beta$ -N-Methylamino-L-alanine in Environmental Samples

J. Beri, K. I. Kirkwood, D. C. Muddiman, et al.

Repligen Corp. now owns the life sciences PAT product portfolio of 908 Devices Inc. Please contact Repligen for further inquiries.

*Anal Bioanal Chem*, Just Accepted Manuscript

DOI: doi.org/10.1007/s00216-018-0930-0

Publication Date (Web): February 17, 2018

Copyright © 2018 Springer-Verlag GmbH Germany, part of Springer Nature

## Contact

Repligen Corporation  
685 Route 202/206  
Bridgewater, NJ, USA 08807  
analytics-support@repligen.com  
(908) 707-1009

## Abstract

We describe a set of new tools for the detection and quantification of  $\beta$ -N-methylamino-L-alanine (BMAA) which includes a novel stable isotope-labeled BMAA standard ( $^{13}\text{C}_3, ^{15}\text{N}_2$ ) and a chip-based capillary electrophoresis mass spectrometry platform for separation and detection. Baseline resolution of BMAA from its potentially confounding structural isomers *N*-2-aminoethylglycine (AEG) and 2,4-diaminobutyric acid (2,4- DAB) is achieved using the chip-based CE-MS system in less than 1 min. Detection and linearity of response are demonstrated across > 3.5 orders of dynamic range using parallel reaction monitoring (PRM). The lower limit of detection and quantification were calculated for BMAA detection at 40 nM (4.8 ng/mL) and 400 nM (48 ng/mL), respectively. Finally, the strategy was applied to detect BMAA in seafood samples purchased at a local market in Raleigh, NC where their harvest location was known. BMAA was detected in a sea scallop sample. Because the BMAA/stable isotope-labeled  $^{13}\text{C}_3, ^{15}\text{N}_2$ -BMAA (SIL-BMAA) ratio in the scallop sample was below the limit of quantification, a semiquantitative analysis of BMAA content was carried out, and BMAA content was estimated to be approximately 820 ng BMAA/1 g of wet scallop tissue. Identification was verified by high mass measurement accuracy of precursor (< 5 ppm) and product ions (< 10 ppm), comigration with SIL-BMAA spike-in standard, and conservation of ion abundance ratios for product ions between BMAA and SIL-BMAA. Interestingly, BMAA was not identified in the free protein fraction but only detected after protein hydrolysis which suggests that BMAA is tightly bound by and/or incorporated into proteins.

[Click here](#) to view the entire article.

