

Design Features of Pre-Packed Disposable, Development to Production Scale Chromatography Columns Confer Reliable Flow and Chromatographic Characteristics

Dana C. Pentia, William J. Wilde, Fletcher L. Malcom, James R. Peyser
Repligen Corporation, Waltham, MA, USA

Summary

Disposable manufacturing has been a developing trend in biopharmaceutical manufacturing for the last two decades. In a 2011 Survey of Biopharmaceutical Manufacturing Capacity and Production¹ the authors note "we are on the verge of the first approvals of mainstream biopharmaceutical products manufactured using single use/disposable systems" and "Single use and disposable chromatography units will be the most important trend in the next five years"².

Disposable technologies have been adopted by biopharmaceutical industry as a mean for faster product changeover, favorable economics, and improved safety. Until now there has not been a broadly applicable solution for disposable chromatography steps.

Repligen's OPUS™ (Open Platform User Specified) pre-packed columns are disposable and can be packed with almost any bioprocessing resin, at internal diameters ranging from 0.5 cm to 60 cm, and column heights from 5 cm and up, offering a broad, scalable solution for the purification of biological products.

Design of the flow distribution permits even flow for all column sizes.


OPUS™ columns characteristics are maintained during various transportation conditions, making them suitable for use at different manufacturing sites.



Design and Materials of Construction

Objective: Design of a simple, clean, robust and functional pre-packed disposable chromatography column

Figure 1: Cross-section of a large scale OPUS™ column

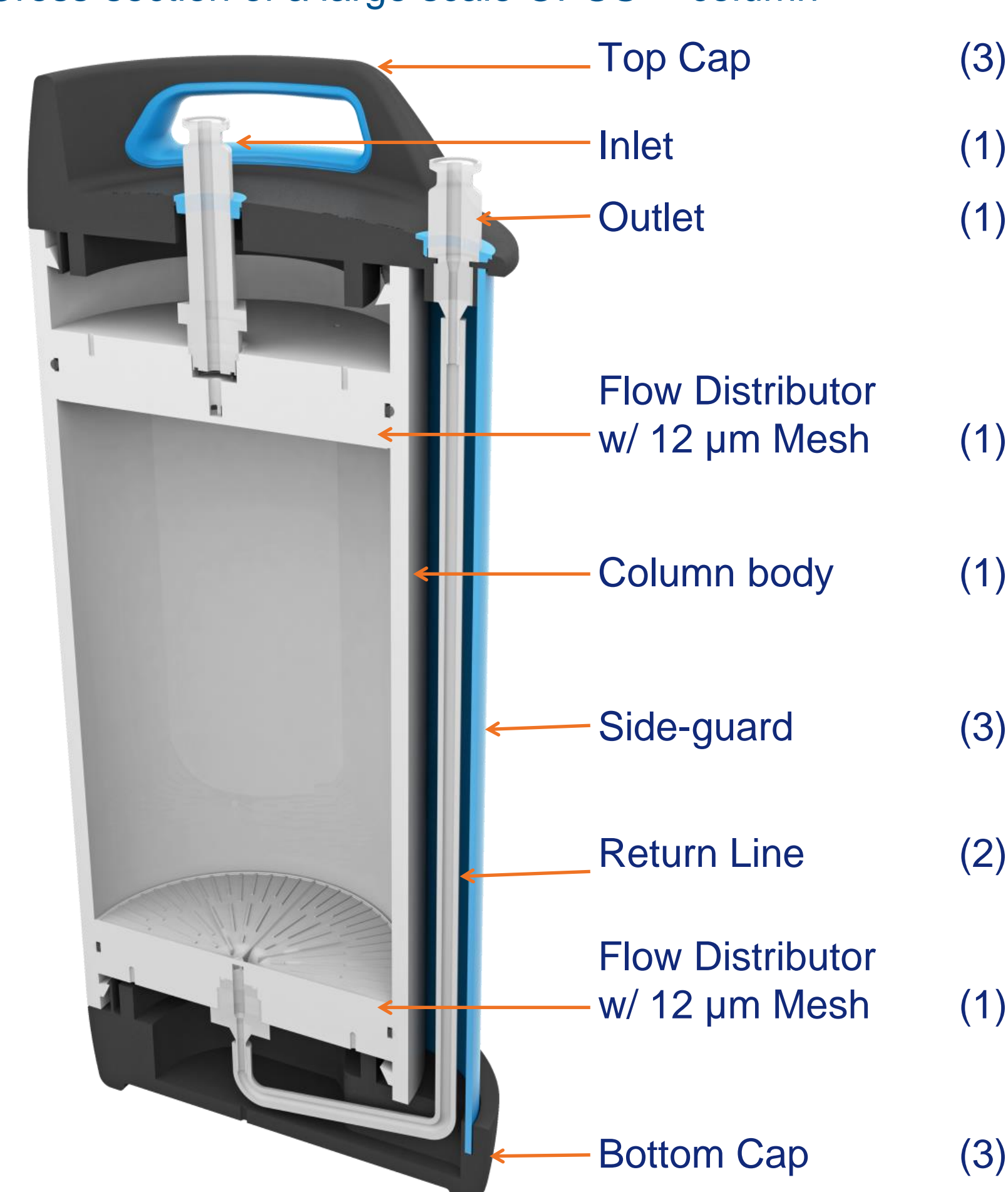

Legend:

Product Contact Material:

- 1 = Polypropylene
- 2 = Platinum Cured Silicone

Non-product Contact Materials:

- 3 = Acrylonitrile Butadiene Styrene (ABS) copolymer

Conclusions:

- Engineered design enables even flow distribution of the mobile phase
- Ease of use is conferred by simplicity of construction
- Regulatory compliance is addressed by the use of USP class VI certified product contact materials

Figure 2: Design features of the flow adaptor

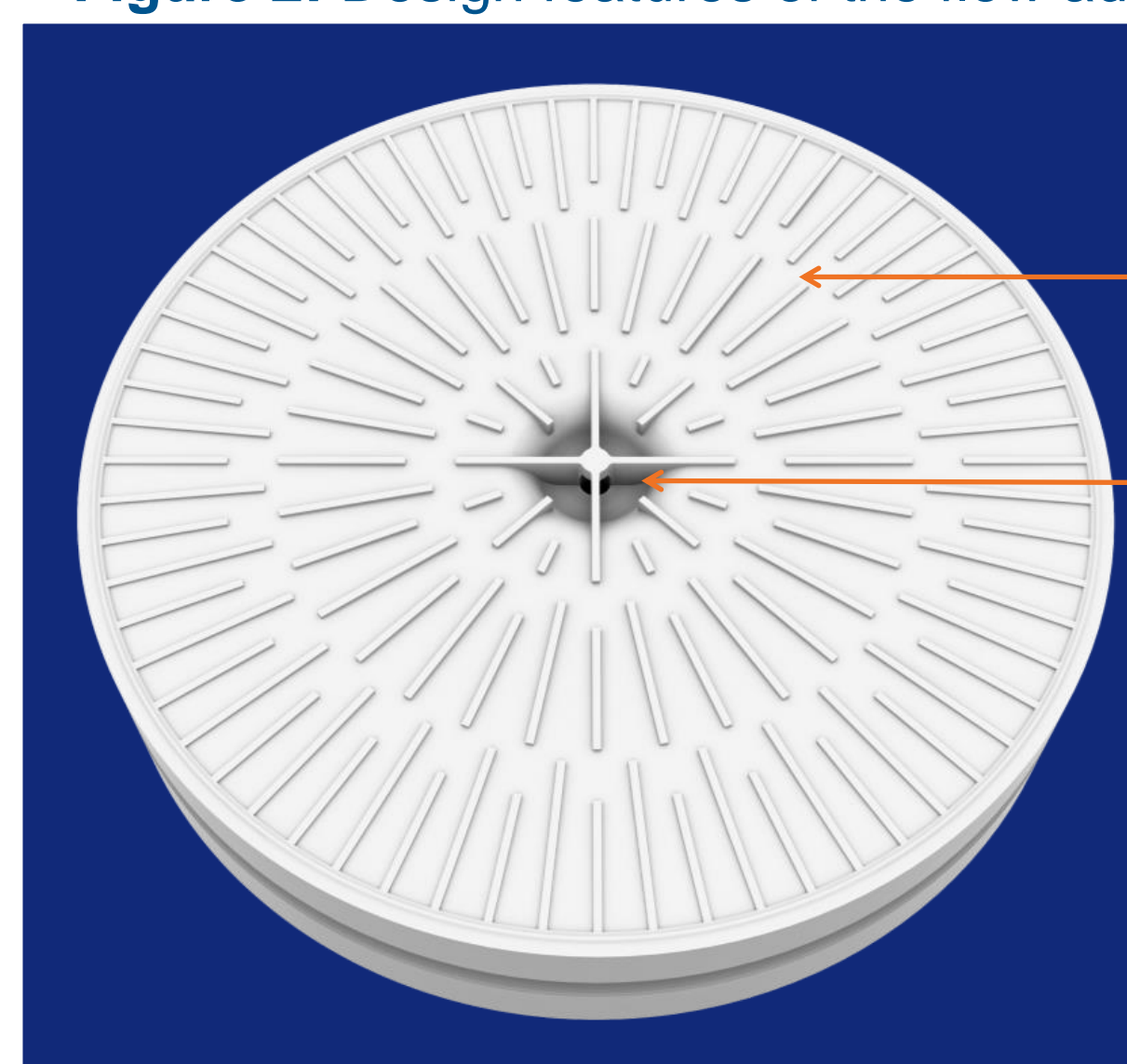


Figure 3: Cross-section of the flow adaptor


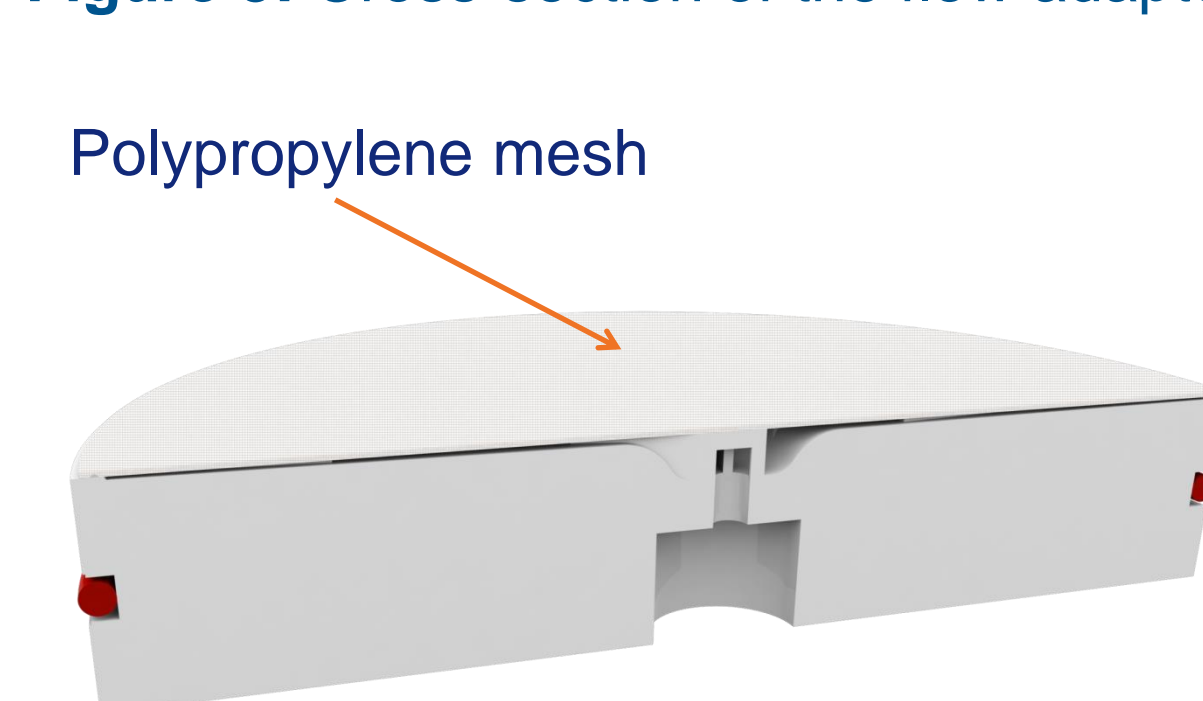


Figure 4: Computational flow mapping

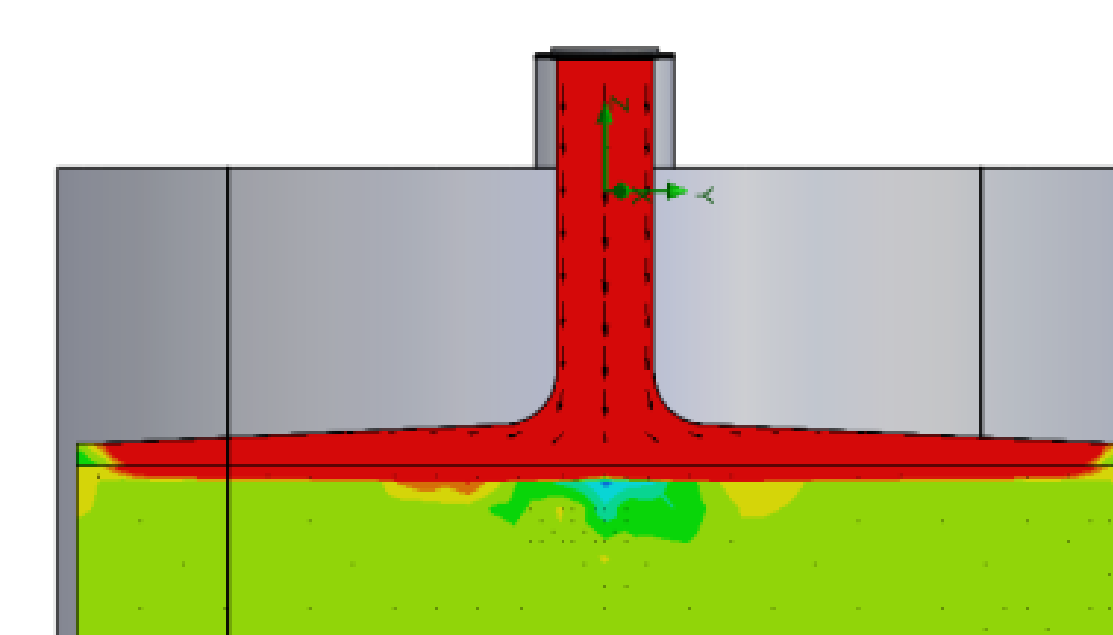

Assessment of Chromatographic Performance

Objective: To experimentally demonstrate flow uniformity of OPUS™ columns through size exclusion performance

Method:

- Determination of plates number and asymmetry for different sizes of OPUS™ columns. Columns are injected with 1% of CV of 2% acetone solution at a linear flow rate of 100 cm/h. Columns are pre-packed with Sepharose® 6FF resin. Columns are all 20 cm height, and IDs of 1.2, 2.5, 8 and 20 cm
- Measuring separation resolution of molecular weight markers on different sizes OPUS™ columns. 1% of CV of 2mg/mL solutions of molecular weight markers injected into the columns at a linear flow rate of 30 cm/h

Acceptance Criteria:

- Column quality attributes pass acceptance test: Plates/m: >1500, Asymmetry: 0.8-1.6
- Resolution of molecular weight markers is consistent for all column scales

Table 1: Column properties for OPUS™ columns pre-packed with Sepharose® 6FF

Column ID	Plates/m @ 100cm/h	Asymmetry @ 100 cm/h
1.2 cm	3010	1.1
2.5 cm	3465	1.2
8 cm	2959	1.2
20 cm	3295	1.2

Conclusion:

- Consistency of separation resolutions and column properties proves flow uniformity
- Chromatographic performance is demonstrated and maintained over a range of OPUS™ column diameters

Figure 5: Chromatogram for separation of molecular weight markers

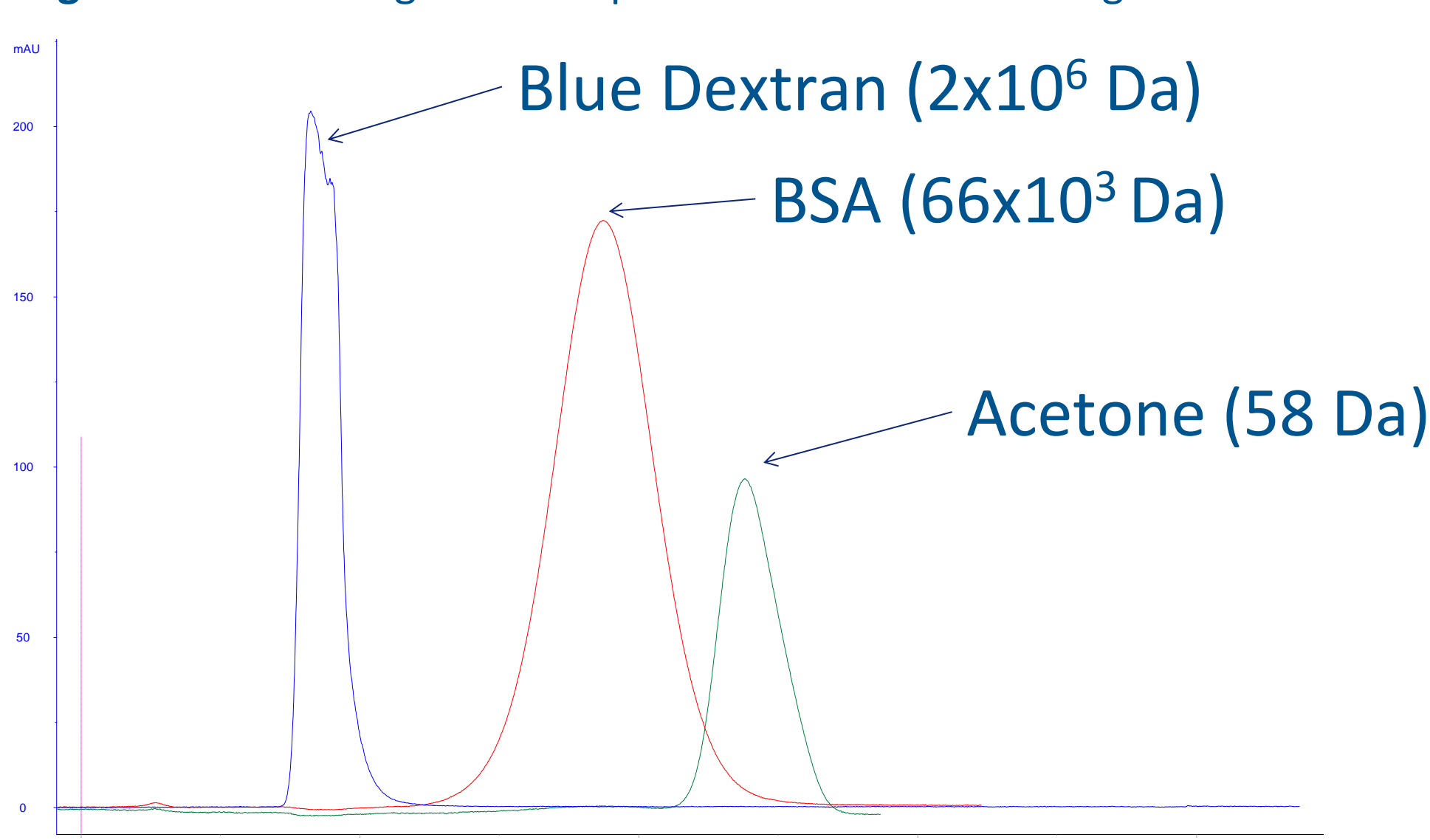


Table 2: Separation resolutions for various sizes OPUS™ columns

Column ID	Resolution Dextran/BSA	Resolution Blue Dextran/Acetone	Resolution BSA/Acetone
1.2 cm	2.2	4.6	0.9
2.5 cm	2.6	6.1	1.1
8 cm	2.3	6.1	1.2
20 cm	2.8	7.1	1.2

Transportation Validation

Objective: To demonstrate that a GMP-ready pre-packed disposable OPUS™ column can be shipped without compromising performance.

Method:

- A 20 cm diameter by 10 cm bed height OPUS™ column was subjected to a rigorous shipping study using the International Safe Transit Association's (ISTA) Procedure 1A
 - Tests included:
 - Fixed displacement vibration test: a total of ~ 15000 vibrational impacts over 64 min
 - Shock-drop test: drop tests on corners, edges, and faces of the packing box from 70 cm (24 inches) height
- A 20x20 cm column packed with SP Sepharose® was shipped 6000 miles by truck and air. The column was assessed for chromatographic efficiency by measuring plates number and asymmetry

Acceptance criteria:

- Intact column upon visual inspection
- Packed bed chromatographic quality attributes (plates number and asymmetry) are preserved

Results:

- Visual Inspection: Minor abrasions caused by vibration testing, some minor damage to the external package cause by drop test on the angles. Column found undamaged and intact. **The package passed ISTA 1A testing**

Figure 6: Column quality testing pre and post shipping; 20x10 cm OPUS™ column


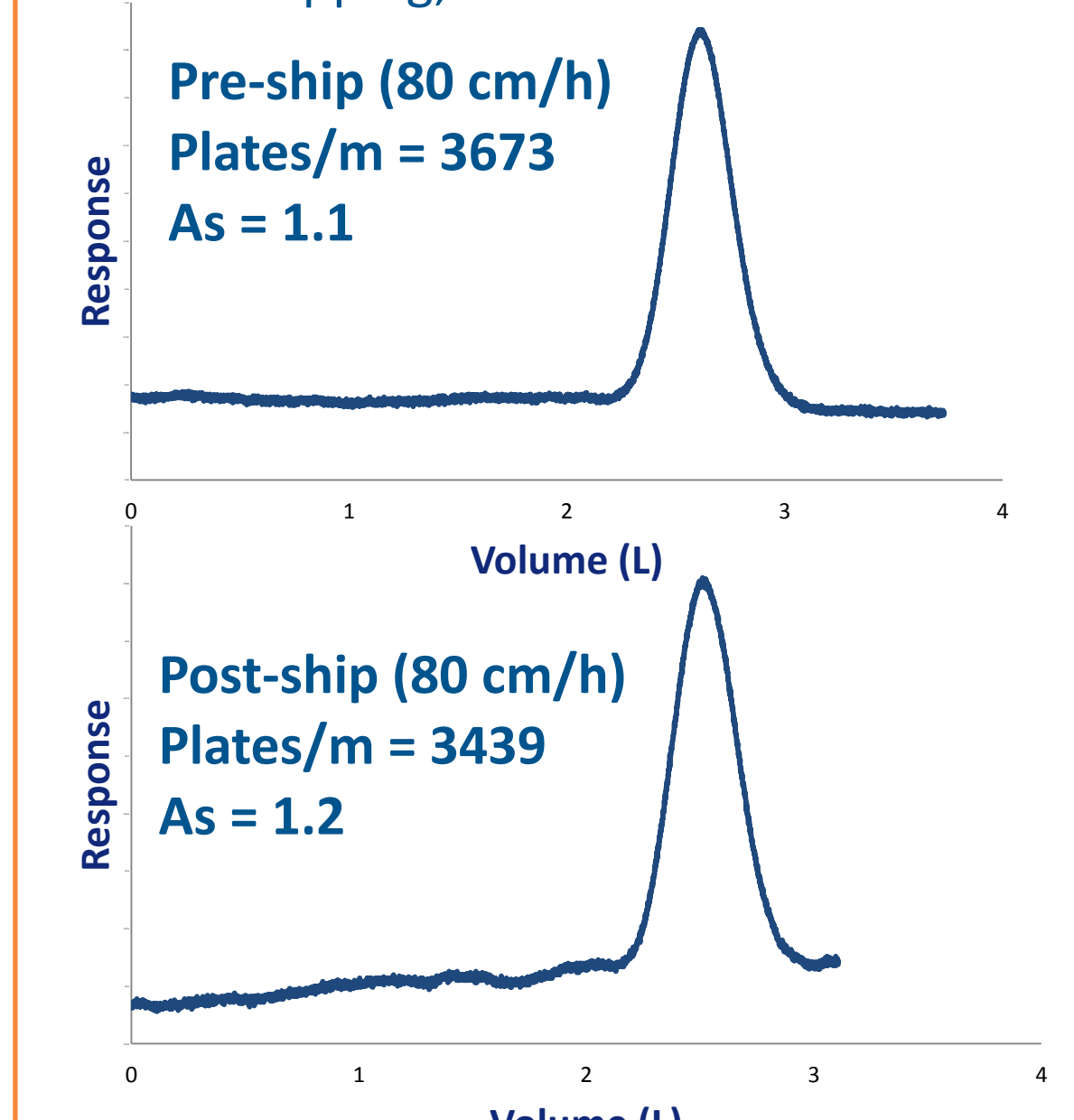


Figure 7: 20x20 cm OPUS™ column quality testing pre and post transportation

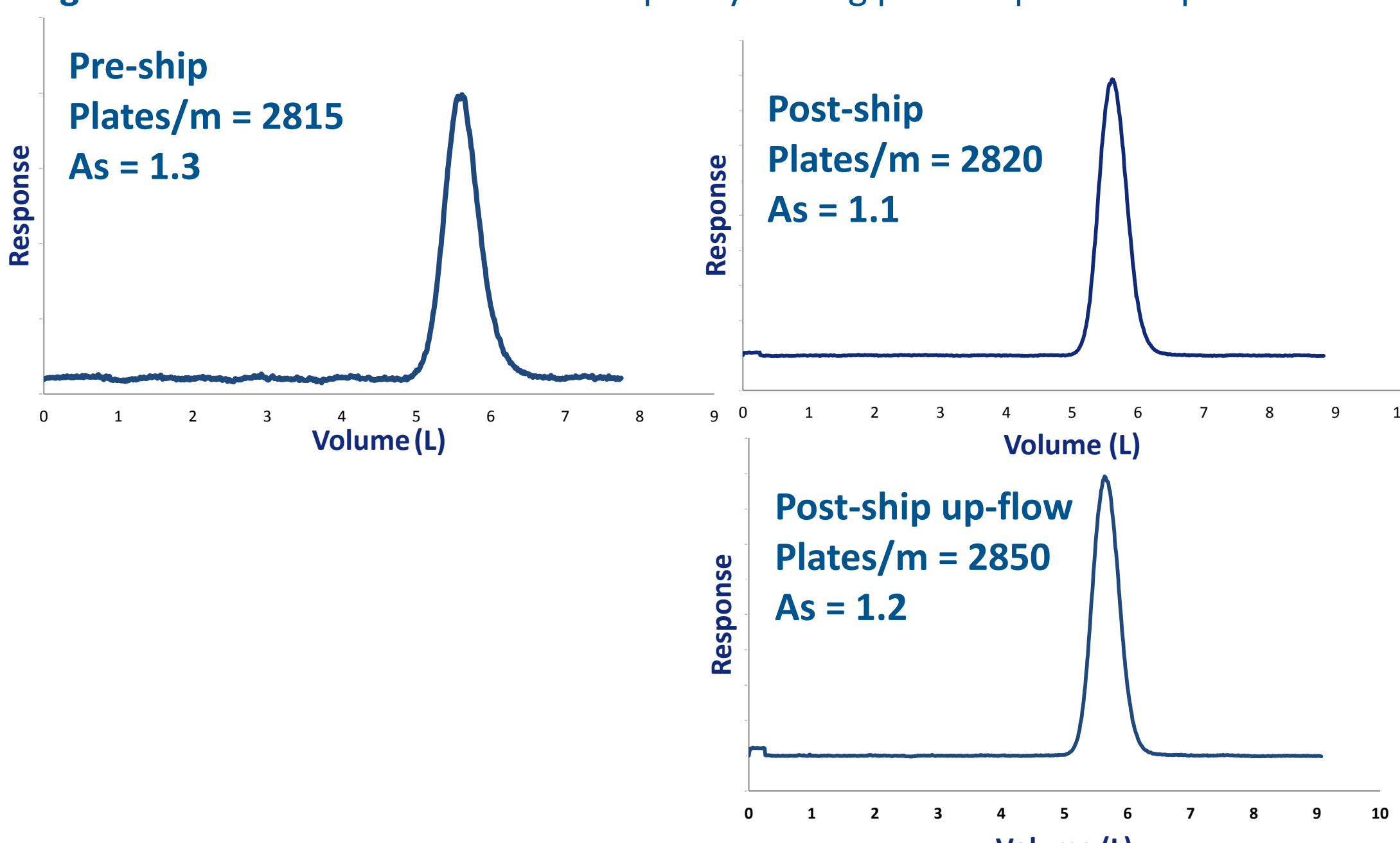

Conclusion:

- Packaging withstood the rigors of a commercial shipping environment
- Chromatographic performance maintained after shipping

Conclusions

- Design and selection of construction materials ensures OPUS™ columns are suitable for manufacturing of biological products
- Engineered design of the flow distributor confers even flow dispersion of the mobile phase
- Outstanding design is confirmed by consistency of size exclusion resolution and column quality attributes for various scales of OPUS™ columns
- Transportability and ruggedness of OPUS™ columns is demonstrated by rigorous tests
- OPUS™ columns offer unparalleled flexibility in regards to column size and choice of chromatographic resin packed, while maintaining the highest standards of quality

Sepharose is a registered trademarks of GEHC
Captiva PriMab is a registered trademark of Repligen Corporation

References:

1. 8th Annual Report and Survey of Biopharmaceutical Manufacturing Capacity and Production
2. Dr. Stefan Schmidt, CSO, ERA Biotech

Company Info

Repligen Bioprocessing
41 Seyon Street Building #1, Suite 100
Waltham, MA 02453
www.repligen.com/bioprocessing